


DRY ICE BUCKET USEAGE TRIALS  
 TRIAL DATE: AUG. 19, 2005

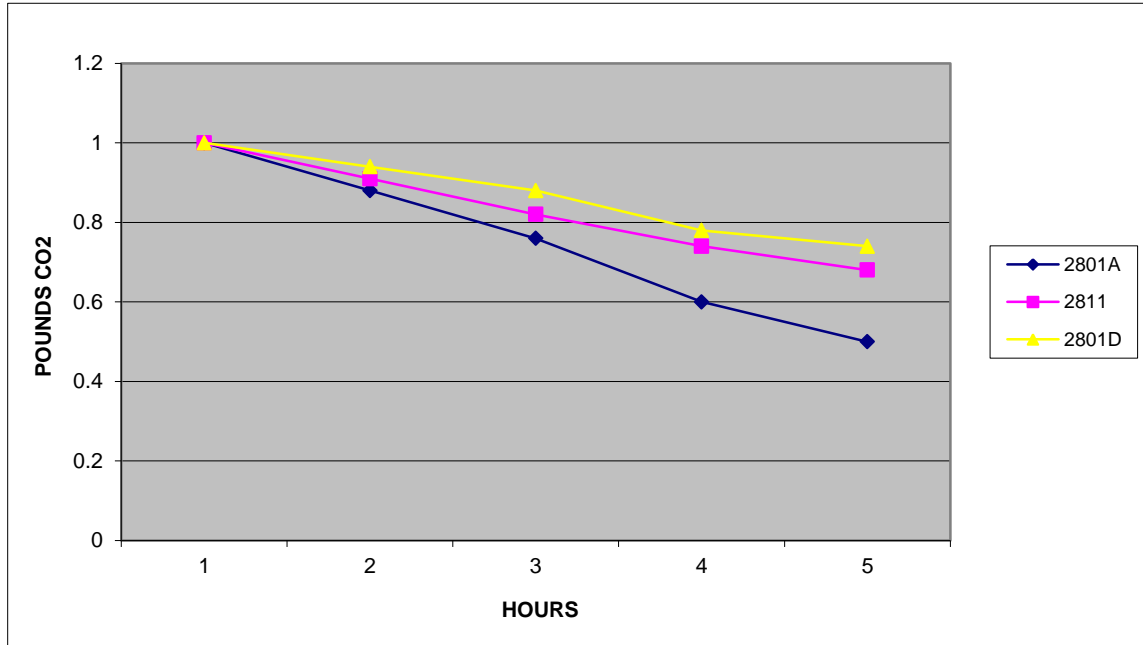

AMBIENT TEMPERATURE 76 DEG.F.

EACH CONTAINER STARTED WITH 1 POUND BLOCK DRY ICE

CONTAINER WEIGHTS IN POUNDS

2801A	0.62	STANDARD 1 GAL.EVS CONTAINER, PLASTAZOTE INSULATION
2811	0.92	1/2 GALLON IGLOO CONTAINER
2801D	1.1	1 GAL. CONTAINER, POLYURETHANE INSULATION

GROSS CONSUMPTION DATA

	1:00 PM START	3:00 PM	4:00 PM	5:00 PM FINISH
2801A	1.62	1.38	1.22	1.12
2811	1.92	1.74	1.66	1.6
2801D	2.1	1.98	1.88	1.84

NET CONSUMPTION DATA

2801A	1	0.88	0.76	0.6	0.5
2811	1	0.91	0.82	0.74	0.68
2801D	1	0.94	0.88	0.78	0.74

CONTAINER #	LBS PER HOUR		HOUR 1	HOUR 2	HOUR 3	HOUR 4
2801A	0.125	2801A	0.12	0.24	0.4	0.5
2811	0.08	2811	0.09	0.18	0.26	0.32
2801D	0.065	2801D		0.12	0.22	0.26