



5/64" Hex Key

Attaching Light Guides To The Eco-20 Light Source

A .280" (7mm) diameter receptacle is provided for attachment of the Fiber optic light guides.

To connect the light guide, simply insert the light guide tip in the receptacle. Insert the 5/64" Hex Key into the set screw (located between the cooling fins, next to the bulb access screw) and tighten. **Do Not Over tighten Set Screw!**

Replacing the Lamp

1. Turn the Eco-20 off and unplug the unit.
2. Remove the thumb screw from the top of the unit and pivot the top cover backward.
3. Remove the 5/64" hex head screw from the bottom of the bulb carrier. Pull out the lamp socket and bulb.


Caution: Do not remove hot lamp with bare fingers.

4. Replace the lamp by sliding it into the socket and pushing the socket back into the bulb carrier. Replace the hex head screw.
5. Pivot the top cover forward and replace the thumb screw.

Lamp Life

The 20 watt lamp has a life of 2000 hours. Allowing unit to cool before moving will extend lamp life.

Replacement bulb - Part #94110

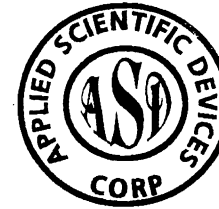

MODEL ECO-20 FIBEROPTIC LIGHT SOURCE

Product Specifications

Kelvin Temperature	3100
Model Number	Eco-20
Lamp Type	94110 (20 watts)
Color	Beige
Height	3"
Width	2 1/2"
Depth	6"
Weight	3 lbs.
Voltage Rating	.2 Amps @ 120 Volts

Parts And Repair Available

From Factory:

- ✕ Replacement lamps
- ✕ Light guides and adapters
- ✕ Repair service

Warrantee

The Eco-20 is guaranteed to be free of manufacturing defects for a period of one year (excludes lamp).

APPLIED SCIENTIFIC DEVICES CORP.
9655 SW VENTURA CT.
TIGARD, OR 97223-9168

503-244-7479
FAX: 503-244-7644